Smart Dilutor W

Instructions for Use

Gravimetric Dilutor

 EN / ES / DE / FR

Doc. 50007662-07

Revision date: April 2026

All Rights Reserved © 2026

CONTENTS

ENGLISH	4
ESPAÑOL	54
DEUTSCH	105
FRANÇAIS	157
IUL Technical Support Contact	209

ENGLISH

Indication of the Introduced Modifications	6
1. Introduction.....	7
1.1. Intended Use of the Smart Dilutor W	7
1.2. General Information	7
2. Safety Information.....	10
2.1. Proper Use	10
2.2. Electrical Safety	10
2.3. Environment	11
2.4. Biological Safety	11
2.5. Chemicals	12
2.6. Maintenance Safety	12
2.7. Waste Disposal	12
2.8. Symbols on the Smart Dilutor W instrument	13
3. General Description.....	14
3.1. Device Overview	14
3.2. Optional Accessories	14
4. Installation	17
4.1. Unpack the Instrument	17
4.2. Site Requirements	17
4.3. Parts Assembly	17
5. Operating Procedures.....	21
5.1. Main screen overview	21
5.2. Placing the bag	24
5.3. Performing a dilution	24
5.4. Performing a dispensing	25
5.5. How to pause or abort	25
5.6. How to create a new program	25
5.7. How to Perform a Calibration	28
5.8. How to Introduce the User ID	29
5.9. How to Establish New Settings	30
6. How to Modify the Stability Function	31
7. External communication and data transfer	32
7.1. Connecting a Barcode Reader	32
7.2. Connecting a PC for data transfer	32
7.3. Connecting a Printer for data transfer	33
7.4. Communication and Output Configuration	34

8.	Maintenance.....	35
8.1.	Cleaning Procedure	35
8.2.	Restoring Factory Settings	35
9.	Troubleshooting.....	37
APPENDIX A	Ordering codes.....	38
APPENDIX B	Technical data.....	41
APPENDIX C	Delivery scope per model.....	44
APPENDIX D	Warranty	51
APPENDIX E	Waste electrical and electronic equipment (WEEE).....	52
APPENDIX F	EU Declaration of Conformity.....	53

Indication of the Introduced Modifications

If it's the first time that you use the Smart Dilutor W, please read the entire document carefully.

If it's not the first time that you use Smart Dilutor W, you can read the indication of the introduced modifications in the table below.

Revision	Description of change
07	Update document format
	Update Section 1 Introduction
	Update Section 1.2.1 Scope of delivery
	Introduce Section 2 Safety Information
	Update Section 3 General Description
	Update Section 4 Installation
	Update Section 5 Operating Procedures
	Introduce Section 7 External communication and data transfer
	Update Section 8 Maintenance
	Update Section 9 Troubleshooting
	Update Appendix A Ordering codes
	Update Appendix B Technical Data
	Update Appendix C Delivery scope per model
	Introduce Appendix D Warranty

1. INTRODUCTION

Thank you for choosing the Smart Dilutor W. We are confident that this instrument will become an integral part of your laboratory.

Before using the Smart Dilutor W, it is essential that you read these Instructions for Use carefully. Following the instructions and safety information will ensure safe operation and maintain the system in a safe condition.

1.1. Intended Use of the Smart Dilutor W

The Smart Dilutor W is a gravimetric diluter designed to perform single-step dilution of samples directly in sterile homogenizer bags and to carry out dispensing operations. It is intended for use with samples that are typically solid, semi-solid, or paste-like, for which gravimetric dilution provides the highest accuracy.

The heart of the Smart Dilutor W is the weighing cell that continuously sends the signal that it generates to the ADC (Analog-to-Digital Converter), then to the microprocessor, which uninterruptedly monitors the weight in the weighing platform and makes the necessary decisions to activate the diluent-pouring pump proportionally. Once the programmed dilution has been reached, the system notifies the user that the process has finished.

NOTE: To obtain optimal performance, the instrument must be kept clean at all times. See the Maintenance Section for more details.

1.2. General Information

1.2.1. Scope of delivery

The Smart Dilutor W comes in several models tailored to meet different workflow needs. Depending on the number of pumps (1 or 2), the type of diluent container (bags or bottles) or whether filters are included (available only for bottle-based models), the package includes specific accessories to ensure optimal setup:

Cat. Number	Description
90007900	SMART DILUTOR W. One pump. For 5 L bottle
90007900BS	SMART DILUTOR W. One pump. For bags
90007901	SMART DILUTOR W. Two pumps. For 5 L bottles
90007901BS	SMART DILUTOR W. Two pumps. For bags
90007901BB	SMART DILUTOR W. Two pumps. For bottles and bags
90007900FI	SMART DILUTOR W. One pump. For 5 L bottle with one filter
90007901FI	SMART DILUTOR W. Two pumps. For bottles with one filter and bags
90007901FII	SMART DILUTOR W. Two pumps. For 5 L bottles with two filters


The delivery of the Smart Dilutor W includes the following items:

For all Smart Dilutor W models:


- A. Smart Dilutor W main unit (Cat. 10007900 or 10007901) with power cord according to destination country).
- B. Arm:
 - **B1**. For 1-pump models (Cat. 90002821).
 - **B2**. For 2-pump models (Cat. 90002822).
- C. Platform and screw for 400 mL bags (Cat. 900012110).
- D. Stainless steel edge of dispensing tube (Cat. 90008331).
- E. Double-sided adhesive tape, 10 units (Cat. 90004955).

For models compatible with **bottles** (Cat. 90007900, Cat. 90007901 & Cat. 90007901BB):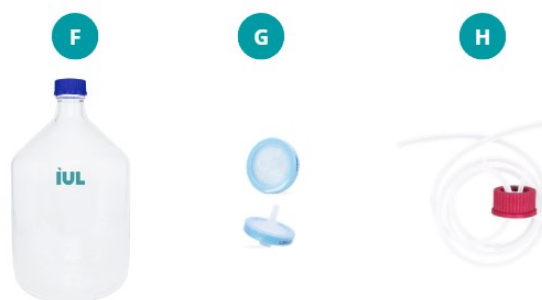

- F. Autoclavable 5 L bottle (GL 45) (Cat. 90008165).
- G. Set of 2 autoclavable 0.2 µm filters (Cat. 90008182).
- H. Tubing set with GL 45 lid (1560 mm) (Cat. 90002825)*.

For models compatible with **bags** (Cat. 90007900BS, Cat. 90007901BS, Cat. 90007901BB):

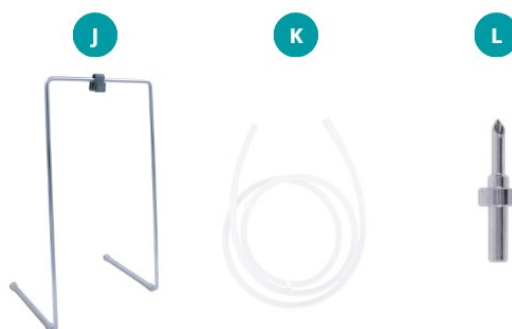- J. Media bag holder (Cat. 90002738).
- K. Silicone tube (1210 mm) (Cat. 90002854)*.
- L. Edge for Septum bags (Cat. 90007528)*.

* Different sizes or lengths are available upon request (see the Optional Accessories section for more details).

For models compatible with **filter(s)** (Cat. 90007900FI, Cat. 90007901FI, Cat. 90007901FII):

- M. Inlet filter tubing (Cat. 900012256)
- N. Outlet filter tubing (Cat. 900014636)
- O. Sterile filter 0.45 μm for bottles (Cat. 90005053)
- P. Filter holder (Cat. 900014621)

Apart from the list of items above, the delivery also comes with the following documents:

- Design qualification (Doc. 50007663)
- Smart Dilutor W Datamatrix Instructions for use (Doc. 50008877)

For a comprehensive list of the scope of delivery per model, including quantities of each item, please refer to Delivery scope per model Appendix.

2. SAFETY INFORMATION

Before using or servicing the Smart Dilutor W, it is essential that you read this user manual carefully. Following the instructions and safety information in this user manual will ensure safe operation and maintain the system in a safe condition.

The following types of safety information appear throughout the Smart Dilutor W Instructions for Use.

WARNING 	The term WARNING is used to inform you about situations that could result in personal injury to you or other people. Details about these circumstances are given in a box like this one.
---	---

CAUTION 	The term CAUTION is used to inform you about situations that could result in damage to the instrument or other equipment. Details about these circumstances are given in a box like this one.
---	--

The advice given in this manual is intended to supplement, not supersede, the normal safety requirements prevailing in the user's country.

2.1. Proper Use

WARNING/ CAUTION 	Risk of personal injury and material damage Improper use of the Smart Dilutor W instrument may cause personal injury or damage to the instrument. The instrument must only be operated by qualified personnel.
--	---

CAUTION 	Damage to the instrument Avoid spilling water or chemicals onto the Smart Dilutor W instrument. Damage caused by water or chemical spillage will void your warranty.
---	---

In case of emergency, switch off the Smart Dilutor W at the power switch and unplug the power cord from the power outlet.

2.2. Electrical Safety

If the operation of the Smart Dilutor W is interrupted in any way (e.g., due to interruption of the power supply or a mechanical error), first switch off the instrument using the power switch, then disconnect the electrical cord from the power outlet. Contact IUL SA Technical Service after such an incident.

WARNING 	Electrical hazard Any interruption of the protective conductor (earth/ground lead) inside or outside the instrument or disconnection of the protective conductor terminal is likely to make the instrument unsafe. Intentional interruption is prohibited. Lethal voltages inside the instrument When the instrument is connected to line power, its terminals may be live. Opening covers or removing parts is likely to expose live parts. In case of replacement or loosening of the mains power cord, this must be replaced <u>only</u> with one of the power cords listed in delivered by IUL SA.
---	---

To ensure safe and satisfactory operation for Smart Dilutor W routine use or after servicing:

- The line power cord must be connected to a line power outlet that has a protective conductor (earth/ground).
- The instrument must not be operated without the cover.
- If liquid has spilled on the instrument, switch it off, disconnect it from the power outlet, and contact IUL SA Technical Services.

If the Smart Dilutor W becomes electrically unsafe, prevent other personnel from operating it, and contact IUL SA Technical Service.

The instrument may be electrically unsafe if:

- The instrument or the power cord appears to be damaged.
- The instrument has been stored under unfavorable conditions for a prolonged period.
- A different mains power cord or fuse is used other than the one provided by IUL SA.

CAUTION 	Risk of explosion or leakage of the internal battery The Smart Dilutor W contains a coin cell battery (CR2032) inside the instrument. The battery must only be replaced by authorized service personnel. Do not attempt to open the instrument to access or replace the battery. Improper replacement may result in damage to the instrument or risk of explosion.
--	--

2.3. Environment

Operating conditions

WARNING 	Explosive atmosphere The Smart Dilutor W is not designed for use in an explosive atmosphere.
---	--

2.4. Biological Safety

Use safe laboratory procedures as outlined in publications such as Biosafety in Microbiological and Biomedical Laboratories, HHS:

https://www.cdc.gov/labs/pdf/SF_19_308133-A_BMBL6_00-BOOK-WEB-final-3.pdf

WARNING 	Samples containing infectious agents Some samples used with the Smart Dilutor W may contain infectious agents. Handle such samples in accordance with the required safety regulations. The responsible person(s) (e.g., laboratory manager) must take the necessary precautions to ensure that the workplace is safe and that the instrument operators are suitably trained and not exposed to hazardous levels of infectious agents, as
---	--

	<p>defined in the applicable Safety Data Sheets (SDSs) or OSHA¹ ACGIH² or COSHH³ documents.</p> <p>Venting of fumes and disposal of wastes must be in accordance with all national, state, and local health and safety regulations and laws.</p>
--	---

2.5. Chemicals

<p>WARNING</p> 	<p>Hazardous chemicals</p> <p>Some chemicals used with the Smart Dilutor W may be hazardous. Always wear safety glasses, gloves, and a lab coat.</p> <p>The responsible person(s) (e.g., laboratory manager) must take the necessary precautions to ensure that the workplace is safe and that the instrument operators are suitably trained and not exposed to hazardous levels of toxic substances (chemical or biological), as defined in the applicable Safety Data Sheets (SDSs) or OSHA, ACGIH or COSHH documents.</p> <p>Venting of fumes and disposal of wastes must be in accordance with all national, state, and local health and safety regulations and laws.</p>
---	--

2.6. Maintenance Safety

<p>CAUTION</p> 	<p>Damage to the instrument</p> <p>Do not use spray bottles containing alcohol or disinfectants to clean the surface of the Smart Dilutor W instrument or its parts. Spray bottles should be used only to clean items that have been removed from the instrument.</p>
---	--

<p>WARNING</p> 	<p>Risk of electric shock</p> <p>Do not open the panels on the instruments.</p> <p>Risk of personal injury and material damage</p> <p>Only perform maintenance that is specifically described in these Instructions for Use.</p>
---	--

2.7. Waste Disposal

Used consumables, such as sample tubes or bags, may contain hazardous chemicals or infectious agents. Such waste must be collected and disposed of properly in accordance with local safety regulations.

For disposal of waste electrical and electronic equipment (WEEE) see the corresponding appendix.

¹ OSHA: Occupational Safety and Health Administration (United States of America).

² ACGIH: American Conference of Government Industrial Hygienists (United States of America).

³ COSHH: Control of Substances Hazardous to Health (United Kingdom).


2.8. Symbols on the Smart Dilutor W instrument


Symbol	Location	Description
	Type plate on the back of the equipment	Waste Electrical and Electronic Equipment (WEEE).
	Type plate on the back of the equipment	CE Mark, Declaration of Conformity, see the corresponding appendix.

3. GENERAL DESCRIPTION

3.1. Device Overview

The device overview of the Smart Dilutor W with two pumps, using diluents in a bottle and a bag (Cat. 90007901BB) is shown below:

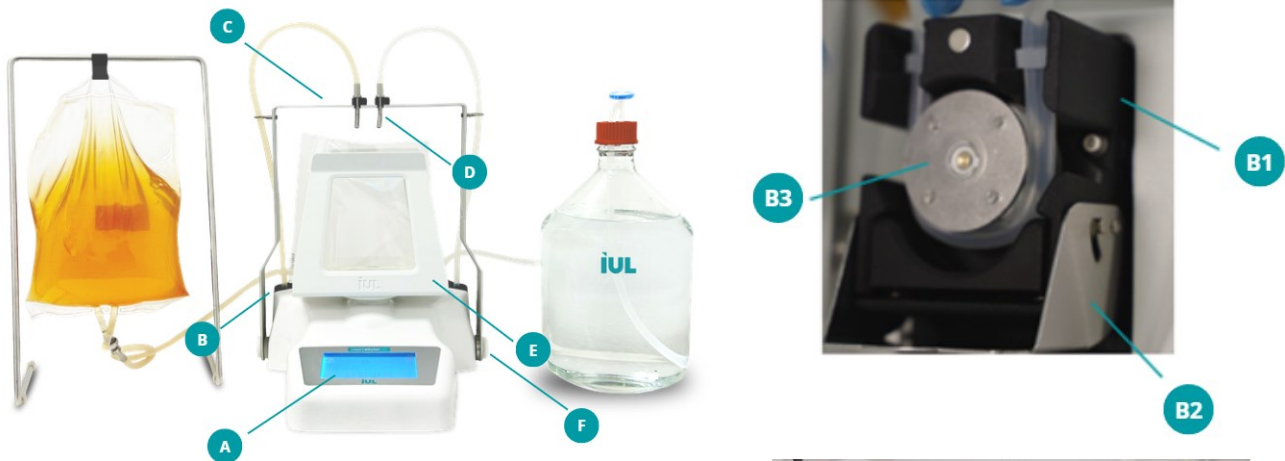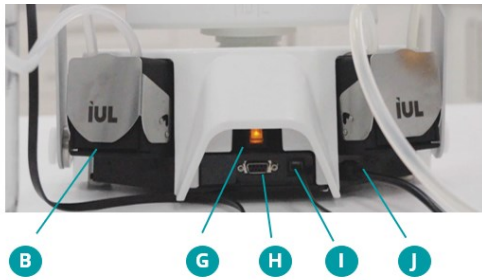- A. Keyboard & Display.
 - B. Pump.
 - B1. Pump head.
 - B2. Pump door.
 - B3. Pump rotor.
 - C. Arm for 2-pump models.
 - D. Stainless steel edge of dispensing tube.
 - E. Platform and screw for 400 mL bags.
 - F. Screw for arm for 2-pump models.
- 
- The bottom view of the device shows callouts G-J: G (Main switch), H (PC and Printer Connection Port (RS-232 DB9M Socket)), I (IUL Barcode reader Connection Port (Special Socket)), and J (Power supply with integrated cable (According to destination country)).
- G. Main switch.
 - H. PC and Printer Connection Port (RS-232 DB9M Socket).
 - I. IUL Barcode reader Connection Port (Special Socket).
 - J. Power supply with integrated cable (According to destination country).

NOTE: The setup of the Smart Dilutor W will vary depending on the model.

For a comprehensive list of the scope of delivery per model, including quantities of each item, please refer to Delivery scope per model appendix.

3.2. Optional Accessories

The Smart Dilutor W is a highly versatile instrument designed to adapt to various laboratory workflows. Some of the components are available in multiple versions to meet specific laboratory requirements.

The tubing sets with GL 45 and silicone tubes provided with the corresponding delivery are also available in extended lengths to offer more flexibility in setup:

- Tubing Set (1645mm) with GL 45 Lid (Cat. 90002897).

- Silicone Tube (1300mm) (Cat. 90002855).
- Silicone Tube (2210mm) (Cat. 90002528).

The edge for septum bags is also available in a larger size for enhanced compatibility with bigger septum bags:

- Edge for Septum bags (big) (Cat. 90008332).

In addition, the Smart Dilutor W also features the following optional accessories to offer enhanced versatility and meet the specific needs of different laboratory workflows.

3.2.1. Multilevel Dilution Arm

The multilevel dilution arm (Cat. 90002806) is an adjustable arm designed to precisely position and align different heights of sample bottles, beakers, or flasks, with the silicone tube dispenser to ensure that the diluent is dispensed accurately and without spills.

Simply unscrew and remove the existing arm located on the right side and install the multilevel dilution arm on the Smart Dilutor W.

3.2.2. Weighing Platform for beakers/flasks

The beakers and flasks platform (Cat. 90002838) is designed to accommodate sample bottles, beakers and flasks, providing a stable and secure setup for liquid dispensing.

Simply unscrew and remove the existing platform (Cat. 900012110) of the Smart Dilutor W and then install and secure the beaker/flask platform.

3.2.3. 80 mL Bag Adapter

The 80 mL bag adapter (Cat. 900014196) is a platform designed to hold 80 mL sample bags in your Smart Dilutor W, ensuring a secure fit and efficient liquid dispensing.

Easily place this bag adapter into the existing platform (Cat. 900012110) of the Smart Dilutor W.

3.2.4. IUL Barcode reader

The IUL barcode reader (Cat. 90005701) streamlines data entry for Smart Dilutor W by enabling easy scanning of Sample ID, Diluent ID, and User ID, eliminating the need for manual input.

Additionally, a support (Cat. 90002836) is available to securely mount the IUL Barcode reader onto the Smart Dilutor W, ensuring stability and convenience during use. See the External communication and data transfer section for details on PC connectivity.

3.2.5. PC Communication cable

The PC communication cable (Cat. 90005518) enables you to export results directly to your computer once the process finishes. See the External communication and data transfer section for details on PC connectivity.

3.2.6. Ticket Dot Matrix Printer

The Ticket Dot Matrix Printer (Cat. 90003069) is an external printer for connecting to the Smart Dilutor W to display and print results on receipts or labels once the process finishes.

NOTE: The RS-232 DB9 cable (Cat. 90005510) required for connecting the Ticket Dot Matrix Printer to the instrument is included with printers purchased directly from IUL. See the External communication and data transfer section for details on PC connectivity.

3.2.7. Ticket Thermal Printer

The Ticket Thermal Printer (Cat. 90003070) is another external printer for connecting to the Smart Dilutor W to display and print results on tickets once the process finishes.

NOTE: The RS-232 DB9 cable (Cat. 90005510) required for connecting the Ticket Thermal Printer to the instrument is included with printers purchased directly from IUL. See the External communication and data transfer section for details on PC connectivity.

3.2.8. Capsule Filter

The Sterile filter 0.45 µm for bottles (Cat. 90005053) is for the sterile filtration of aqueous solutions (e.g., buffers, diluents). It features a 0.45 µm pore size with a glass microfiber pre-filter to improve particulate loading and extend filter life, followed by a nylon membrane for bacteria sterile filtration. The filter reduces the need for frequent preparation of autoclaved diluents, thereby simplifying daily operation. The filter is held in place using the filter holder (Cat. 900014621).

4. INSTALLATION

4.1. Unpack the Instrument

The packaging will depend on the accessories and model of Smart Dilutor W purchased.

The packaging of the Smart Dilutor W can be stored for reuse.

4.2. Site Requirements


Place the Smart Dilutor W instrument on a solid, vibration-free surface. Ideally, a firmly secured bench top, near an earthed/grounded electrical outlet. While operating, make sure that the instrument does not touch other objects. Locate it away from vibration-sensitive instruments. Ensure that there are no special difficulties operating the main switch located at the rear part of the instrument.

CAUTION 	Treat the instrument like a precision balance due to its weighing cell.
---	---

4.3. Parts Assembly

Before switching on the instrument, assemble all parts according to your specific model.

For instance, the following drawing shows the assembly of the Smart Dilutor W with two pumps and using diluents in a bag and in a bottle (Cat. 90007901BB).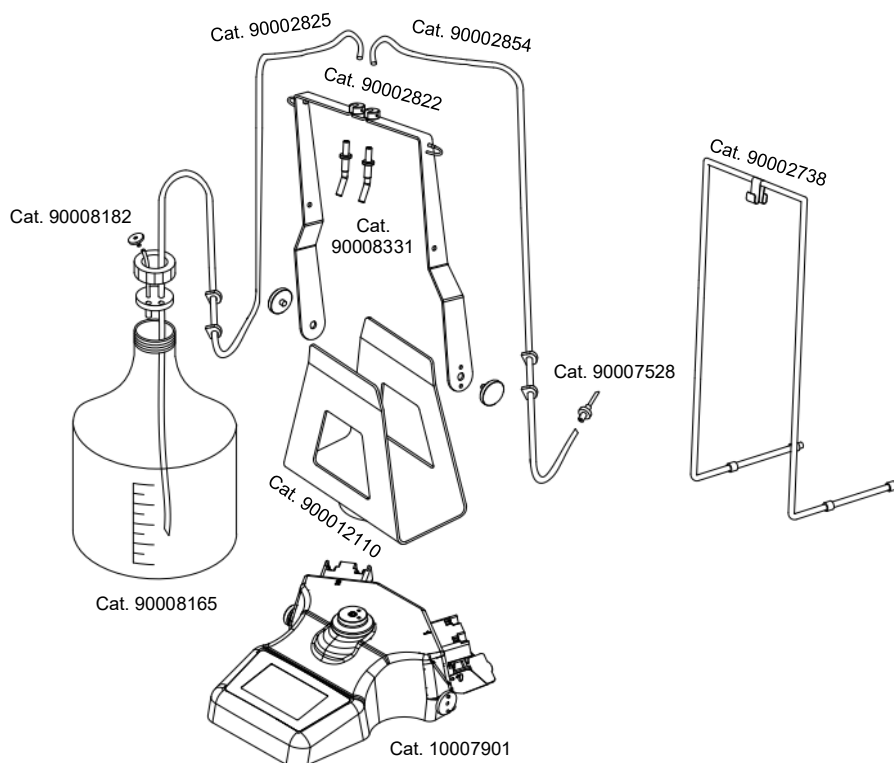

Below is the step-by-step instructions to assemble the instrument:


1. Position the platform (Cat. 900012110) on the weighing cell, hold it in place, and carefully secure it using the inox screw, tighten until fully seated.

NOTE: Ensure it aligns properly with the designated notch. (See the troubleshooting section)


2. Attach the appropriate arm for your model: Cat. 90002821 for 1- models or Cat. 90002822 for 2-pump models.

NOTE: Ensure the arm can move freely back and forth.

3. Open the pump door and carefully place the tube inside. Ensure that the tube clamp (Part B in the photo below) is placed correctly, i.e., above the part encircled in blue in the photo.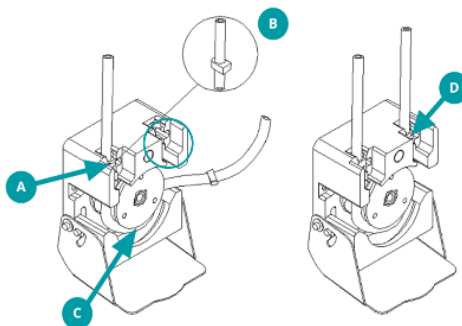

4. Insert the tube into the pump such that the end of the tube coming from the diluent should be placed in the **IN** hole.

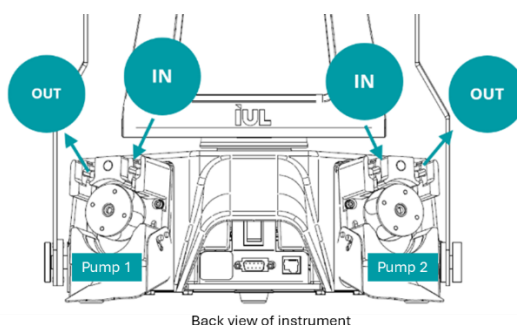Back view of instrument

IN: indicates the hole where the tube must be inserted.

OUT: shows the part where the tube must exit.

NOTE: Pump 1 rotates clockwise. Pump 2 rotates counterclockwise.

NOTE: It's important to place the tube in the intended direction; otherwise, the pump will not be able to move the diluent from the diluent bottle or bag into the sample bag.

5. Position the tube below the rotor.

CAUTION

Avoid twisting the tube 180° to prevent blockages that could obstruct the liquid flow.

6. Guide it out through the opposite hole (**OUT**), securing its corresponding tube clamp in its designated slot.
7. Close the peristaltic head.

NOTE: The tube will be stretched tight, and the pump door can be properly closed. If the tube is not placed correctly, usually due to the incorrect placement of the tube clamp, the tube will be slack, and the pump door cannot be closed completely.

8. Connect the diluent source, either a bottle or a bag, to the tubing system.

- *If working with the diluent in bottles:*

- A. Screw the GL 45 lid of the tubing set (Cat. 90002825) onto an autoclavable bottle (Cat. 90008165).

- B. Attach the 0.2 µm autoclavable filter (Cat. 90008182).

NOTE: The filter prevents contamination by filtering the air entering the bottle.

- *If working with the diluent in bags:*

- A. Place the media bag in the designated media bag holder (Cat. 90002738).

- B. Insert the edge for septum bags (Cat. 90007528) into the silicone tube (Cat. 90002854).

- C. Attach the edge for septum bags (Cat. 90007528) to the media bag.

9. Insert the stainless-steel edge (Cat. 90008331) into the opposite end of the tube and place it into its slot on the arm.

NOTE: Its shape directs flow against the bag wall, minimizing splashes.

10. Plug in the power cord, ensuring the arm is pushed backwards.

11. Turn on the main switch. The instrument will display its firmware version and perform a quick auto-test before reaching the main screen.

4.3.1. Filter Assembly

Follow these steps if you are working with a filter (Cat. 90007900FI, Cat. 90007901FII, Cat. 90007901FI).

Initial setup:

1. Screw the GL 45 lid of the Inlet filter tubing (Cat. 900012256) onto an autoclavable bottle (Cat. 90008165).

2. Connect and insert the open end of the Inlet filter tubing (Cat. 900012256) to the half-stepped barb inlet of the filter.

3. Connect one end of the Outlet filter tubing (Cat. 900014636) to the half-stepped barb outlet of the filter.

4. Attach and insert the stainless-steel edge (Cat. 90008331) to the opposite end of the Outlet filter tubing (Cat. 900014636)

NOTE: There is an arrow on the filter housing that indicates the correct liquid flow direction (toward the outlet) inside the filter. Ensure tubing follows this direction.

5. Follow the steps in Section 4.3 on how to position the Inlet filter tubing (Cat. 900012256) inside the pump.

6. Secure the filter in place using the filter holder (Cat. 900014621). Turn the filter upside down so that the air vent is positioned at the bottom.

Air removal procedure:

1. Instead of placing the filter outlet tubing into a bag, direct it into a suitable container.

2. Do **not** place the container on the Smart Dilutor W platform.

3. Set up a dispensing procedure.

4. Pull the arm forward to start dispensing.

NOTE: Because the container is not placed on the Smart Dilutor W, dispensing will continue indefinitely unless manually stopped.

5. Allow dispensing to continue until the outlet flow is smooth (i.e., no bursts of air bubbles).
6. Push the arm backward to stop dispensing.
7. Turn the filter upright so that the air vent is at the top. A large air bubble may be visible inside the filter.
8. Pull the arm forward to start dispensing again.
9. Slightly open the air vent and tilt the filter toward you so that the air bubble moves directly beneath the vent.

NOTE: Liquid will flow out of the air vent. This is normal and necessary to minimize air remaining inside the filter.

10. Close the air vent.
11. Push the arm backwards to stop dispensing.
12. Turn the filter upside down again, positioning the air vent at the bottom.

System verification:

1. Position the filter outlet tubing appropriately.
2. Place a bag on the Smart Dilutor W.
3. Test the system by performing a standard dispensing procedure.

NOTE: For the Smart Dilutor W with filter, the delivered volume (in dispensing or dilution mode) may show a maximum deviation of ± 0.5 g. This deviation is caused by residual air bubbles within the filter. As additional dispensing or dilution cycles are performed, the deviation will decrease as remaining air bubbles are eliminated.

4. The system is now ready for use.

Between workdays, it is essential to check for air bubbles inside the filter. If the deviation in delivered volume exceeds 0.5 g, repeat the Air Removal Procedure to ensure that air inside the filter is minimized.

When the system is not in use—especially for extended periods—apply the clamp to prevent backflow of liquid, which could introduce air into the filter.

5. OPERATING PROCEDURES

5.1. Main screen overview

Depending on the process (dilution or dispensing), the main screen will be the following:

The screen is divided into three horizontal regions:

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5.1.1. Top region

The top region displays information related to the weight measurements.

Display:	Name:	Function:
	Bag symbol	Displayed when performing a dilution process.
	Bottle symbol	Displayed when performing a dispensing process.
	Tare button	To perform a tare.
	Shaking weight icon	Indicates that the instrument is tolerant to vibrations after the stability function has been set to a value other than the default "0".
	Current weight display	Shows the current weight.

5.1.2. Middle region

The middle region displays information related to the sample.

DISPENSING

DILUTION

Display:	Name:	Function:
<i>Dispensing</i>		
	Filled bottle symbol	Represents the sample.
	Sample ID	Can be introduced manually or by using the IUL Barcode reader connected to the Smart Dilutor W. NOTE: Alphanumeric information can be accepted.
	Dispensing weight	Shows the weight to be dispensed.
	User ID	Shows the user ID.
<i>Dilution</i>		
	Shrimp icon	Represents the sample.
	Sample ID	Can be introduced manually or by using the IUL Barcode reader connected to the Smart Dilutor W. NOTE: Alphanumeric information can be accepted.
	Current sample weight taken	Shows the final stabilized sample weight when the arm is backwards.
	Dilution factor	Shows the dilution factor used.
	User ID	Shows the user ID.

5.1.3. Bottom region


The bottom region displays information related to the workflow settings.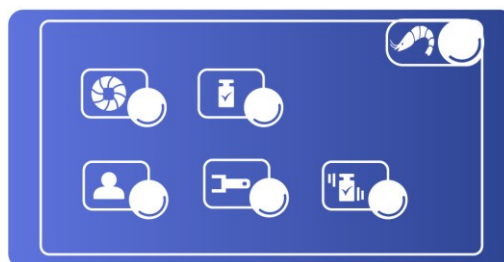

Display:	Name:	Function:
	Arm icon and triangle	A flashing triangle icon is shown briefly indicating the position of the arm:
		backwards (upwards triangle) or
		forward (downwards triangle).
	Pump icon	Shows the pump number currently in use (1 or 2, if applicable).
	Diluent ID	This ID is optional.
	Utilities button	Opens the utilities dialog menu

5.1.4. Utilities dialog menu

To set more configurable parameters, the Utilities button can be press and the following the Utilities dialog menu will be displayed:

Display:	Name:	Function:
	Pump icon	Establish a new program (see the 'How to create a new program' section for detailed instructions).
	User icon	Establish the User ID (see the 'How to introduce the user ID' section for detailed instructions).
	Calibration weight icon	Perform calibration (see the 'How to perform a calibration section for detailed instructions).
	Spanner icon	Establish the User ID (see the 'How to establish new settings' Section for detailed instructions).
	Shaking weight icon	Establish a new stability function (see the 'How to modify the stability function' section for detailed instructions).
	Shrimp icon and Flask icon	Both icons allow you to return to the previous screen. The Shrimp icon is displayed on the top-right side of the screen during the dilution process, while the Flask icon appears during the dispensing process.

5.2. Placing the bag

1. Align the bag mouth with the upper edge of the platform.

NOTE: The upper part of the platform has double-sided adhesive tapes (Cat. 90004966) to secure the bag in place.

2. Gently bring the sides of the platform together until they grip the mouth of the bag.
3. Release the platform and ensure the bag remains open and ready for filling.

5.3. Performing a dilution

1. The bag symbol () briefly flashes to indicate the procedure is underway.
2. Press the pump button () to choose the pump (1 and/or 2, if applicable).
3. Press the tare button () to cancel out the bag weight. It flashes when the instrument asks for a tare action.

NOTE: During the taring process, ---.- is displayed on the current weight display ().

4. Carefully add the sample to the bag.

NOTE: The Shrimp icon () flashes while the unit waits for the sample.

NOTE: The grams (g) character flashes if the weight measurement isn't stable.

5. Gently pull the arm forward.
6. Observe the pump dispensing the diluent until reaching the target weight; the current weight displays on the screen.

NOTE: The pump slows down as it nears the target weight.

7. A short beep can be heard when the target weight is reached.

NOTE: If the ticket printer is correctly connected, a ticket will be printed.

Push the arm backwards to prepare for a new dilution or keep the arm forward if another dilution is ready to be done.

5.4. Performing a dispensing

1. The Bottle symbol () briefly flashes to indicate that the procedure is underway.
2. Press the pump button () to choose the pump (1 and/or 2), if applicable.
3. Place the container (bag or bottle) where the diluent must be dispensed.
4. Press the Tare button () to cancel out the container weight.

NOTE: During the taring process, “---.” is displayed on the current weight display ().

NOTE: The Filled-bottle icon () flashes while the unit awaits the diluent.

5. Gently pull the arm forward.
6. Observe the pump dispensing the diluent until reaching the target weight; the current weight displays on the screen.
NOTE: The pump slows as it nears the target weight.
7. A short beep can be heard when the target weight is reached.
NOTE: If the ticket printer is connected, a ticket will be printed.
8. Push the arm backwards to prepare for a new dispensing.

5.5. How to pause or abort

Pausing or aborting can be performed in two ways:

- A. Pushing the arm tube backwards allows you to pause the operation (dilution or dispensing process). Pulling the arm forward again will resume the operation until reaching the target weight.
NOTE: When pushing the arm backwards, do so gradually until the motor stops. This gradual movement helps prevent spillage from the bag. Once the motor stops, push the arm backwards completely.
- B. Pressing the pause button will also pause the operation. Simply press it again to resume the process.

5.6. How to create a new program

After selecting the Pump icon on the Utilities dialog menu, the following options will be available:


Choose between a dispensing (Flask icon) or a dilution (Shrimp icon):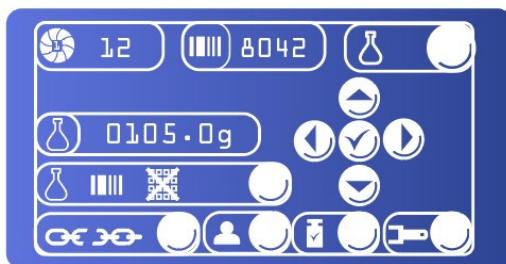


- **Dispensing:** Click the Flask icon () to set up a dispensing program.
- **Dilution:** Click the Shrimp icon () to set up a dilution program.

5.6.1. New dispensing

After selecting the Flask icon, the following dispensing screen will be displayed:

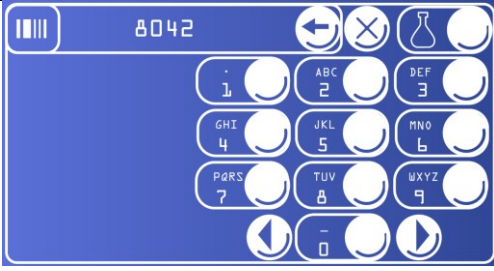
Display:	Name:	Function:
	Pump icon	<p>If applicable, choose the pump to be used for the dispensing process by navigating with the left and right arrows.</p> <p>NOTE: Optionally, once the pump is selected, you can decide whether to continue with dilution or switch to the dispensing process by using the up and down arrows. If dispensing is selected, the screen will automatically switch to the dilution settings menu.</p> <p>The Smart Dilutor W stores the last-used program per pump and per mode:</p> <ul style="list-style-type: none"> - 2 programs with a 1-pump unit (1 for dispensing + 1 for dilution). - 4 programs with a 2-pump unit (2 for dispensing + 2 for dilution).
	Diluent ID	<p>Set the diluent ID.</p> <p>NOTE: Can be introduced manually or by using the IUL Barcode connected to the unit.</p>
	Flask icon with weight	Specify the weight to be dispensed.
	Arrows and check button	Used to navigate and confirm displayed values
	Input method button	<p>Select input method (keyboard or IUL Barcode reader). If the barcode reader is installed, it will light up when an operation starts, and the barcode will be stored as the Sample ID.</p> <p>NOTE: Only the last 5 digits will display on the screen.</p> <p>NOTE: The default sample ID is a number that increases automatically with each operation.</p> <p>NOTE: The keyboard icon is tagged by default; press its near-right button to remove the tag and switch to keyboard input for each new operation. The main screen will display up to 6 digits.</p>

	
 	<p>Chain icon</p> <p>Allows the used pump 1 to work with pump 2 either together (closed chain) or separately (broken chain). This option is displayed when two pumps are installed.</p> <p>NOTE: When the closed chain is selected, the pump number that will work with the current pump is shown.</p>

5.6.2. New dilution

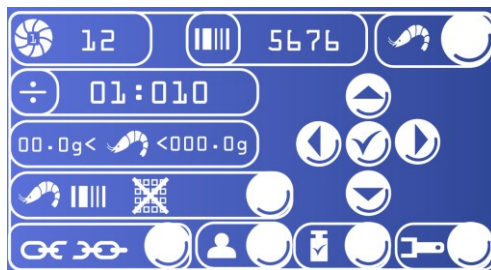After selecting the Shrimp icon () , the following dilution screen will appear:

Display:	Name:	Function:
	Pump icon	<p>If applicable, choose the pump to be used for the dispensing process by navigating with the left and right arrows.</p> <p>NOTE: Optionally, once the pump is selected, you can decide whether to continue with dilution or switch to the dispensing process by using the up and down arrows. If dispensing is selected, the screen will automatically switch to the dilution settings menu.</p> <p>So, the Smart Dilutor W stores the last-used program per pump and per mode:</p> <ul style="list-style-type: none"> - 2 programs with a 1-pump unit (1 for dispensing + 1 for dilution). - 4 programs with a 2-pump unit (2 for dispensing + 2 for dilution).
	Diluent ID	<p>Set the diluent ID.</p> <p>NOTE: Can be introduced manually or by using the IUL Barcode connected to the unit.</p>

	Dilution factor	Specify the dilution factor.
	Limits of the sample weight	Define the sample weight limits. If the weight sample is not within these limits, the dilution process will not be performed.
	Arrows and check button	Used to navigate and confirm displayed values
	Input method button	<p>Select input method (keyboard or IUL Barcode reader). If the barcode reader is installed, it will light up when an operation starts, and the barcode will be stored as the Sample ID.</p> <p>NOTE: Only the last 5 digits will display on the screen.</p> <p>NOTE: The default sample ID is a number that increases automatically with each operation.</p> <p>NOTE: The keyboard icon is tagged by default; press its near-right button to remove the tag and switch to keyboard input for each new operation. The main screen will display up to 6 digits.</p>
	Chain icon	<p>Allows the used pump 1 to work with pump 2 either together (closed chain) or separately (broken chain). This option is displayed when two pumps are installed.</p> <p>NOTE: When the closed chain is selected, the pump number that will work with the current pump is shown.</p>

5.7. How to Perform a Calibration

<p>CAUTION</p>	<p>Calibration should be conducted on a stable surface devoid of vibrations, with the stability function set to "0" (default value) and with the bag holder installed.</p>
-----------------------	--

After selecting the Calibration weight icon () on the Utilities dialog menu, the following options will be available:

Display:	Name:	Function:
	Tare button	Used to zero the weight.
	Calibration weight icon	Used to initiate the calibration process.
	Calibration weight arrow keys	Use the arrows to fine-tune the calibration weight. The current measured weight is displayed in the lower box.

Follow these steps to perform a calibration

- Tare button check** (): If it is flashing, press it to perform a tare. Upon completion, the Calibration weight icon () will begin flashing.
- Positioning the calibration weight:** Place the calibration weight at the center of the platform.
- Adjustment** (if necessary): Use the arrows to fine-tune the calibration weight (upper box -) according to the calibration weight used. The current measured weight is displayed in the lower box ()
- Initiating Calibration:** Press the Calibration weight icon () to initiate the calibration process.
NOTE: If the weight appears incorrect, or if calibration cannot be completed successfully, please contact IUL for assistance.
NOTE: If the Ticket Dot Matrix Printer (Cat. 90003069) or Ticket Thermal Printer (Cat. 90003070) is connected, a ticket will be printed.

5.8. How to Introduce the User ID

After selecting the User icon () on the Utilities dialog menu, use the arrows and confirmation button to change the values:

- Manual Entry:** Modify the User ID using the keyboard.


- **Barcode Reader Entry:** Use the IUL Barcode reader to input the User ID.

NOTE: Alphanumeric information is accepted.

NOTE: The main screen will display only the last 8 digits.
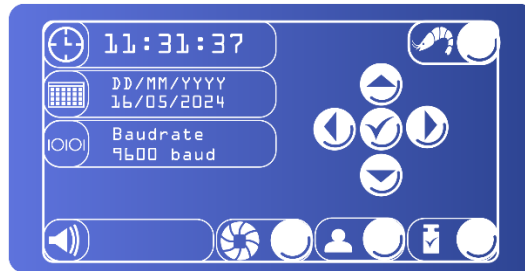
5.9. How to Establish New Settings

After selecting the Spanner icon () on the Utilities dialog menu, the following options will be accessible:

Display:	Name:	Function:
 11:31:37	Time setting	Configure the current time.
 DD/MM/YYYY 16/05/2024	Date setting	Configure the current date.
 Baudrate 9600 baud	Printer/PC communication parameters	Adjust communication settings (see Section 7.3 for detailed instructions).
	Sound Settings	Enable or disable the end-operation buzzer.

6. HOW TO MODIFY THE STABILITY FUNCTION

After selecting the Shaking weight icon () on the Utilities dialog menu, the following screen will appear:

This function enables you to adjust the tolerance of surface vibrations where the Smart Dilutor W is positioned.

NOTE: Placing the instrument on a mechanical vibrating surface may result in a significant delay in reaching the operation endpoint.

Select stability level: use the arrows to choose the desired stability level:

- **“0” (default):** suitable for stable surfaces with minimal vibration.
- **“4”:** highest level, suitable for surfaces experiencing strong vibrations.

<p>CAUTION</p> 	<p>Choosing higher stability levels on stable surfaces may reduce dilution time but compromise accuracy.</p>
---	--

7. EXTERNAL COMMUNICATION AND DATA TRANSFER

The following peripherals can be connected to the back of the Smart Dilutor W:

- IUL Barcode Reader (Cat. 90005701)
- PC
- Ticket Dot Matrix Printer (Cat. 90005510) and Ticket Thermal Printer (Cat. 90003070).

7.1. Connecting a Barcode Reader

Only IUL barcode reader (Cat. 90005701) is compatible with the Smart Dilutor W.

Connect the barcode reader to the dedicated terminal located on the instrument.

To configure the IUL barcode reader, see Barcode Reader Instructions For Use (Doc. 50005710).

When the instrument prompts for a Sample ID, Diluent ID, or User ID on the display, the scanned code is automatically entered into the corresponding field.

7.2. Connecting a PC for data transfer

Connect the PC communication cable to the RS-232 DB9M socket located on the back of the Smart Dilutor W and to your computer's USB port.

Once connected, the communication parameters must be configured in the Smart Dilutor W menu as described in the 'Communication and Output Configuration' section.


The results are generated in real time and are not stored in the instrument, so the PC must remain connected during operation.

Every operation records the following information: Software version (e.g. IUL Smart Dilutor 2.20.), Date and time, Operator ID, Sample ID, Sample Weight, Dilution Factor, Diluent ID, Total Weight, Final Accuracy and Last Calibration.

To transfer data from Smart Dilutor W, you can use the Windows software Hyperterminal. If it's unavailable in your Windows version, consider the free Termiter software as an alternative.


Other terminal programs can be used to set up the same communication parameters.

1. Open the Terminal Software: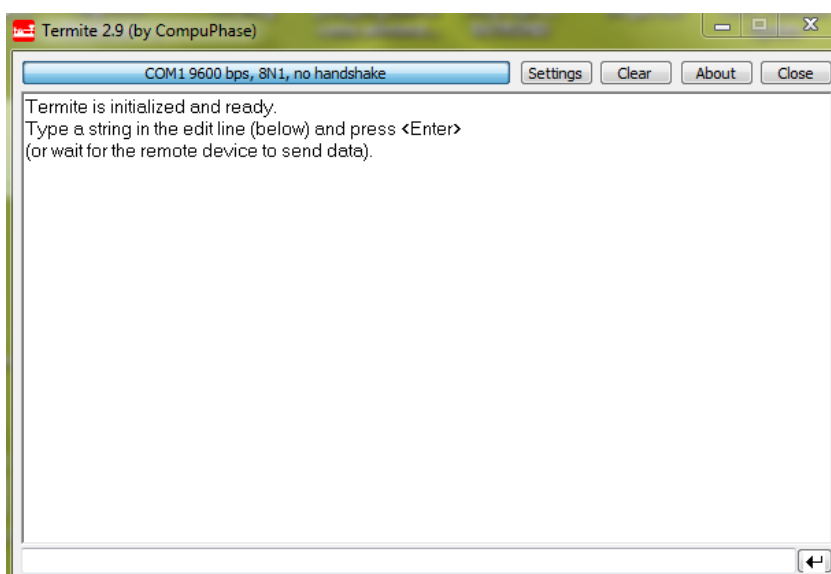


2. On the Terminal Software select the Computer COM port where the Smart Dilutor W is connected and set the following communication parameters:

3. Click on settings
 - Baud rate: 9600
 - Data bits: 8
 - Parity: Even
 - Stop bits: 1
 - Flow control: RTS/CTS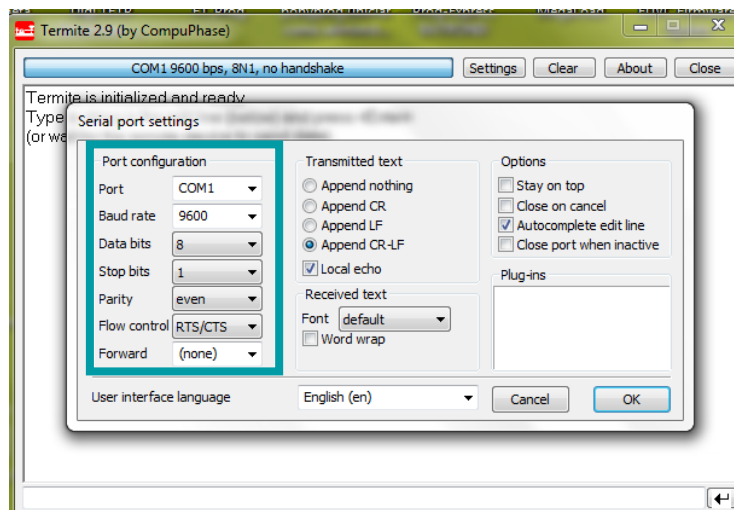

4. Click "OK" on the Terminal Software. The instrument sends the data in this format, and it is displayed as follows: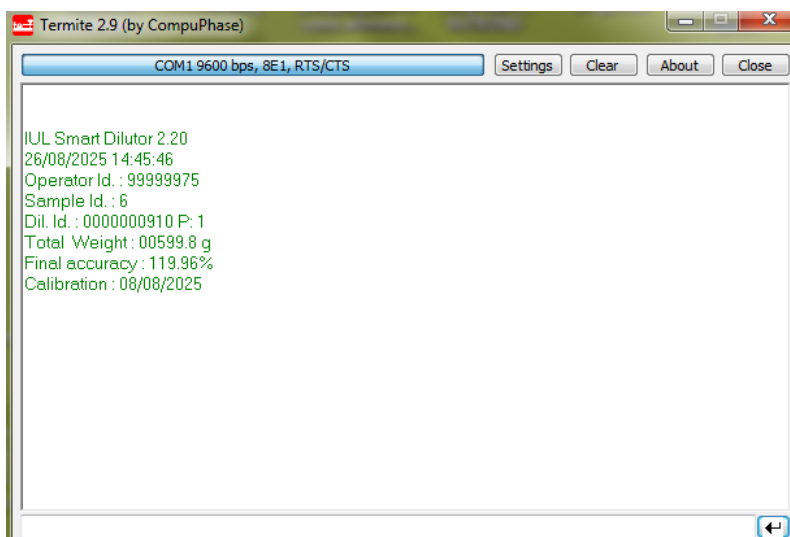

This information can be copied and pasted into a third-party software such as Microsoft Excel or Microsoft Word.

7.3. Connecting a Printer for data transfer


Printers must be connected using the serial cable supplied with the printer (Cat. 90005510).

Once connected, the communication parameters must be configured in the Smart Dilutor W menu as described in the following section.

The results are generated in real time and are not stored in the instrument; they cannot be retrieved retrospectively, so the printer must remain connected during operation.

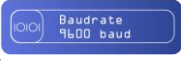
Every operation records the following information: Software version (e.g. IUL Smart Dilutor 2.20.), Date and time, Operator ID, Sample ID, Sample Weight, Dilution Factor, Diluent ID, Total Weight, Final Accuracy and Last Calibration.

Example of printed ticket for Ticket Dot Matrix Printer:

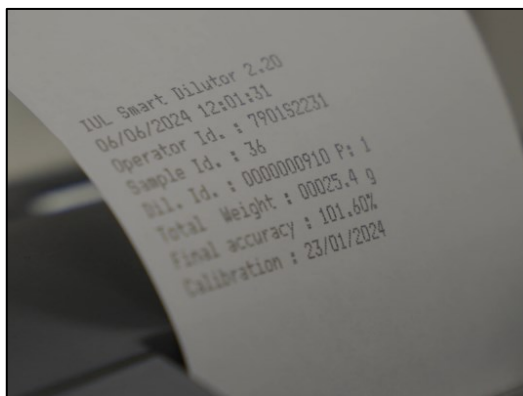7.4. Communication and Output Configuration

The following steps show how to configure the required communication parameters of the Smart Dilutor W to ensure proper communication with the connected PC or printer.

1. Press the Spanner icon () on the Utilities dialog menu to adjust communication settings.
2. Go to Printer or PC communication parameters (, Baudrate 9600 baud).
3. In 'RS Interface', select 'enable' option:
 - **Enable:** Requests different parameters to be set:
 - **Baudrate settings:** Set 9600 baud.
 - **Data bits:** Set 8 bits.
 - **Parity:** Set even.
 - **Stop bits:** Set 1 or 2 bits.
 - **Flow Ctrl:** Set RTS/CTS

NOTE: Select "Disabled" option if no RS Interface is connected.
4. If required 'Com Format', it requests different parameters to be set:
 - Data Format: Set the structure of the data. Choose 'CSV' or 'Printer'.
 - If CSV is set, it requests a Separator:
 - Separator: To set the character to be used when separating data from archives. Set 'Semicolon', 'Space', 'Tab', 'User Def.' or 'Comma'.

NOTE: User Def. is required to be written down.

8. MAINTENANCE

8.1. Cleaning Procedure

Before cleaning, disconnect the Smart Dilutor W from the power supply.

Clean the Smart Dilutor W main unit, arm, platform and screw for 400 mL bags, media bag holder, edge for septum bags and stainless-steel edge of the dispensing tube using a cloth moistened with water and soap. Dry the bag platform screw and the platform support nut before replacing the platform.

To disinfect these components, wipe the external surfaces with a cloth moistened with deionized water and 70% ethanol or isopropyl alcohol (except for the platform for 400 mL bags).

Clean the platform for 400 mL bags using water with 3% bleach.

<p>WARNING</p> 	<p>A special cleaning of the bag holder screw and the support nut shall be performed once per month using an IPA solution to prevent potential damage to both components. Ensure that the inner surface of the support nut is thoroughly cleaned and completely dry before reassembly. An IPA swab may be used to clean and dry the inner area of the support nut on the bag holder.</p> 
---	--

Clean the screen using a soft brush, hot water, and soap applying careful movements to avoid scratching the surface.

Clean the tube with water and soap, then perform one dispensing cycle with clean water to rinse the internal path.

The following components can be autoclaved:

- Silicone tube
- Tubing set with GL 45 lid
- Edge for septum bags
- Stainless-steel edge of the dispensing tube
- Autoclavable 5 L bottle (GL 45)
- Set of 2 autoclavable 0.2 µm filters

8.2. Restoring Factory Settings

Error or malfunctions may be caused by wrong parameter values.

Restore the factory settings to recover the standard value for every parameter and let the instrument be in a known “must run” state.

Beware that the PC and printer settings are also reset.

After restoring factory settings when a printer is used, the communication parameters must be reset.

To restore factory settings:

1. Turn off the instrument.

2. Turn it on again, and when the splash screen appears, keep the upper left zone pressed until the default parameters dialog appears:
3. Use the up and down arrows to select the desired value (Yes/No).
4. Press the Shrimp icon to return to the main menu.

9. TROUBLESHOOTING

Issue	Cause	Solution
Both pumps work at the same time unintentionally	Enchained pumps	Select the unchained pumps  . See the 'How to create a new program' section.
When calibrating, the weight is different in each side of the bag holder	The calibrated weight is not placed in the center of the bag holder	Repeat the calibration with the calibrated weight in the center of the bag holder
When calibrating, the calibration weight is not detected	The bag holder is not placed	Place the bag holder, and then perform the calibration
Error Pump 1 or Error Pump 2	The motor in pump 1 or 2 is stuck	Clean the pump head as described in the 'Maintenance' section.
Error Pump 1 or Error Pump 2	The motor in pump 1 or 2 is stuck	Check that the tube is placed correctly, in the right direction, and the tube is tight. Refer to instructions indicated in the 'Parts assembly' section.
The liquid in the tube goes back to the bottle or media bag	The pump header doesn't press well the tube	Check the tube for damages
The liquid in the tube goes back to the bottle or media bag	The pump header doesn't press the tube enough	Check that the tube is placed correctly, in the right direction, and the tube is tight. Refer to instructions indicated in the 'Parts assembly' section.
The bag holder is not straight or cannot be screwed to the support.	The bag holder is not well placed.	Check the support where the bag holder is placed as well as the bottom pin of the bag holder. Must be like that: 

APPENDIX A ORDERING CODES

Spare parts

Product	Contents	Cat.
Arm for 1-pump models	1 unit	90002821
Arm for 2-pump models	1 unit	90002822
Double-sided adhesive tape	10 units	90004955
Stainless steel edge of dispensing tube	1 unit	90008331
Platform and screw for 400 mL bags	1 unit	900012110
Inox screw + washer	1 unit	900012157

Spare parts (for Bottles)

Product	Contents	Cat.
Tubing set (1560 mm) with GL 45 lid	1 unit	90002825
Autoclavable 5 L bottle (GL 45)	1 unit	90008165
Set of 2 autoclavable filters 0.2 µm	1 unit	90008182

Spare parts (for Bags)

Product	Contents	Cat.
Media bag holder	1 unit	90002738
Silicone tube (1210 mm)	1 unit	90002854
Edge for Septum bags	1 unit	90007528

Spare parts (for Bottles with filters)

Product	Contents	Cat.
Sterile filter 0.45 µm for bottles	1 unit	90005053
Filter holder	1 unit	900014621
Inlet filter tubing	1 unit	900012256
Outlet filter tubing	1 unit	900014636
Filter kit for bottles	1 unit	90002898
<i>(1 unit each: Sterile filter 0.45 µm for bottles (90005053), Filter holder (900014621), Inlet filter tubing (900012256), and Outlet filter tubing (900014636))</i>		

Optional accessories

Product	Contents	Cat.
Multilevel dilution arm	1 unit	90002806
Support for IUL Barcode reader	1 unit	90002836
Weighing platform for beakers/flasks	1 unit	90002838
80 mL bag adapter	1 unit	900014196
Ticket Dot Matrix Printer (Cat. 90005510 included)	1 unit	90003069
Ticket Thermal Printer (Cat. 90005510 included)	1 unit	90003070
Printer cable	1 unit	90005510
IUL Barcode reader	1 unit	90005701

Optional accessories (for Bottles)

Product	Contents	Cat.
Tubing set (1645mm) with GL 45 lid	1 unit	90002897

Optional accessories (for Bags)

Product	Contents	Cat.
Silicone tube (1300mm)	1 unit	90002855
Silicone tube (2210mm)	1 unit	90002528
Edge for Septum bags (big)	1 unit	90008332

APPENDIX B TECHNICAL DATA

IUL SA reserves the right to change specifications at any time.

WEIGHING SPECIFICATIONS

Weight resolution	0.1 g (tenth of a gram)
Weight accuracy	≥ 99 % (for weight ≥ 5 g)
Calibration weight	100 - 2000 g, user configurable (100 g steps)
Surface vibration compensation levels	0 - 4 (5 different levels)

DILUTION SPECIFICATIONS

Dilution ratio	XX (Sample) / XXX (Sample + Diluent)
Maximum total weight	2400 g
Sample weight resolution	0.01 g (hundredth of a gram)
Minimum sample weight	1 g
Maximum sample weight	Depends on the dilution ratio (within the maximum weight capacity).
Limits of the sample weight	User configurable, from 00.0 g to 999.9 g The dilution process will not be performed if the sample weight is outside these limits
Dilution accuracy	≥ 99% (for dilution weight ≥ 10 g)
Dilution deviation*	± 0.5 g (independent of delivered weight)
Delivery speed	~ 16 mL/s

*for filter models ONLY

DISPENSING SPECIFICATIONS

Dispensing range	5 – 1000 g
Dispensing accuracy	≥ 99% (for weight ≥ 10 g)

Dispensing deviation* ± 0.5 g (independent of delivered weight)

Delivery speed ~ 16 mL/s

*for filter models ONLY

NOTE: Accuracy specifications comply with the requirements of ISO 7218:2007/Amd.1:2013 and ISO 6887-1:1999.

GENERAL FEATURES

Dimensions (LxWxH)* 34 x 32 x 37 cm / 13.4 x 12.6 x 14.6 in

Weight* 6.5 kg / 14.3 lb

Connectivity PC and Printer Port (RS-232 DB9M Socket).

IUL Barcode Reader Port (special Socket).

Touch screen interface 5"

Program memory Last-used program per pump and mode (dispensing/dilution) — up to 4 programs (2- pump unit) / up to 2 programs (1-pump unit).

End-of-operation alarm Yes (enable or disable)

** Total dimensions/weight of Smart Dilutor W 1-pump model main unit - Cat. 10007900, Arm for 1-pump models - Cat. 9000282, Platform and screw for 400 mL bags - Cat. 900012110*

OPERATING CONDITIONS

Mains 100 - 240 ACV 50/60 Hz

Maximum power 40 W

RTC Backup Battery Coin Type Lithium Manganese Dioxide Battery

Model CR2032 3 Volt, 220mAh

Overvoltage category II

Air temperature 15 °C – 32 °C / 41°F – 89 °F

Relative humidity 15 – 75 % (non-condensing)

Altitude Up to 2000 m / 6500 ft

Place of operation	For indoor use only
Pollution level	2

TRANSPORTATION CONDITIONS

Air temperature	-25 °C – 60 °C / -13 °F – 140 °F In the manufacturer's packaging
Relative humidity	Maximum of 75 % (non-condensing)

STORAGE CONDITIONS

Air temperature	5 °C – 40 °C / 41 °F – 104 °F In the manufacturer's packaging
Relative humidity	Maximum of 85% (non-condensing)

PACKAGE DIMENSIONS AND WEIGHT

Cat.	Dimensions (LxWxH)	Weight
90007900	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	10.5 kg / 23.1 lb
90007900BS	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	9.5 kg / 20.9 lb
90007901	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	14.5 kg / 32.0 lb
90007901BS	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	10.0 kg / 22.0 lb
90007901BB	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	12.0 kg / 26.5 lb
90007900FI	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	10.6 kg / 23.3 lb
90007901FII	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	14.8 kg / 32.6 lb
90007901FI	60 x 40 x 40 cm / 23.6 x 15.7 x 15.7 in	12.1 kg / 26.7 lb

APPENDIX C DELIVERY SCOPE PER MODEL

The description of the delivery scope for each Smart Dilutor W model is detailed here. The delivery scope is divided into two tables, a table for the models with 1 pump and the second table for the models with 2 pumps. For each model, the quantity delivered per item is indicated.

Smart dilutor W with 1 pump


One Pump		For Bottle	For Bag	For Bottle with 1 filter
		90007900	90007900BS	90007900FI
Item	Image			
Smart Dilutor W 1-pump model main unit (Cat. 10007900)		1 unit	1 unit	1 unit
Arm for 1-pump models (Cat. 90002821)		1 unit	1 unit	1 unit
Platform and screw for 400 mL bags (Cat. 900012110)		1 unit	1 unit	1 unit
Stainless steel edge of dispensing tube (Cat. 90008331)		1 unit	1 unit	1 unit

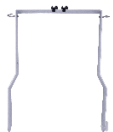Double-sided adhesive tape (Cat. 90004955)		10 units	10 units	10 units
Autoclavable 5 L bottle (GL 45) (Cat. 90008165)		1 unit	N/A	1 unit
Tubing set (1560 mm) with GL 45 lid* (Cat. 90002825)		1 unit	N/A	N/A
Set of 2 autoclavable 0.2 µm filters (Cat. 90008182)		1 unit	N/A	1 unit
Media bag holder (Cat. 90002738)		N/A	1 unit	N/A
Silicone tube (1210 mm) (Cat. 90002854)*		N/A	1 unit	N/A
Edge for Septum bags (Cat. 90007528)*		N/A	1 unit	N/A
Sterile filter 0.45 µm for bottles (Cat. 90005053)		N/A	N/A	1 unit

Filter Holder (Cat. 900014621)		N/A	N/A	1 unit
Inlet filter tubing (Cat. 900012256)		N/A	N/A	1 unit
Outlet filter tubing (Cat. 900014636)		N/A	N/A	1 unit

* Different sizes or lengths are available upon request (see 'Optional Accessories' Section).

Smart Dilutor W models with two pumps

Two Pumps		For Bottles	For Bags	For Bottle + Bag	For Bottles with 2 Filters	For Bottle with 1 Filter + Bag
		Cat. 90007901	Cat. 90007901BS	Cat. 90007901BB	Cat. 90007901FII	Cat. 90007901FI
Item	Image					
Smart Dilutor W 2-pump model main unit (Cat. 10007901)		1 unit	1 unit	1 unit	1 unit	1 unit
Arm for 2-pump models (Cat. 90002822)		1 unit	1 unit	1 unit	1 unit	1 unit
Platform and screw for 400 mL bags (Cat. 900012110)		1 unit	1 unit	1 unit	1 unit	1 unit
Stainless steel edge of dispensing tube (Cat. 90008331)		2 units	2 units	2 units	2 units	2 units

Double-sided adhesive tape (Cat. 90004955)		10 units	10 units	10 units	10 units	10 units
Autoclavable 5 L bottle (GL 45) (Cat. 90008165)		2 units	N/A	1 unit	2 units	1 unit
Tubing set (1560 mm) with GL 45 lid* (Cat. 90002825)		2 units	N/A	1 unit	2 units	1 unit
Set of 2 autoclavable 0.2 µm filters (Cat. 90008182)		2 units	N/A	1 unit	2 units	1 unit
Media bag holder (Cat. 90002738)		N/A	1 unit	1 unit	N/A	1 unit
Silicone tube (1210 mm) (Cat. 90002854)*		N/A	2 units	1 unit	N/A	1 unit
Edge for Septum bags (Cat. 90007528)*		N/A	2 units	1 unit	N/A	1 unit
Sterile filter 0.45 µm for bottles (Cat. 90005053)		N/A	N/A	N/A	2 units	1 unit

Filter Holder (Cat. 900014621)		N/A	N/A	N/A	2 units	1 unit
Inlet filter tubing (Cat. 900012256)		N/A	N/A	N/A	2 units	1 unit
Outlet filter tubing (Cat. 900014636)		N/A	N/A	N/A	2 units	1 unit

* Different sizes or lengths are available upon request (see 'Optional Accessories' Section).

APPENDIX D WARRANTY

The Smart Dilutor W instrument has a 12-month warranty period which starts when the instrument is purchased from IUL SA.

The warranty is void when misuse of the equipment can be proved. Damage or manufacturing faults caused by impacts, chemical or corrosive products, liquids, dampness, or other external factors, such as radiation, fire, or inadequate transport are not included.

In addition, the warranty will not apply if the equipment has been handled, repaired, or modified by unqualified or specifically designated personnel.


As part of this warranty, all shipping costs to the selected technical service are the responsibility of the Buyer.

APPENDIX E WASTE ELECTRICAL AND ELECTRONIC EQUIPMENT (WEEE)

This section provides information about the disposal of waste electrical and electronic equipment by users.

The crossed-out wheeled bin symbol (see below) indicates that this product must not be disposed of with other waste; it must be taken to an approved treatment facility or to a designated collection point for recycling, according to local laws and regulations.

The separate collection and recycling of waste electronic equipment at the time of disposal helps to conserve natural resources and ensures that the product is recycled in a manner that protects human health and the environment.

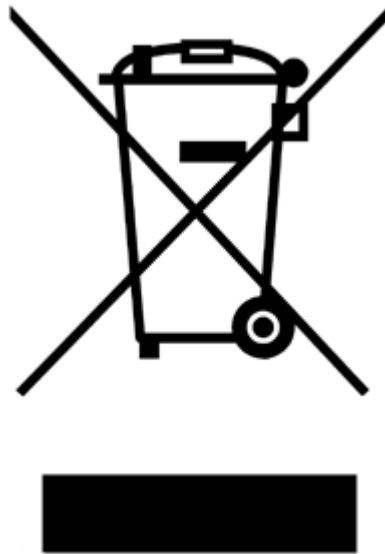APPENDIX F EU DECLARATION OF CONFORMITY

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Cat. 90007901

Conforms the Council Directives:

- Directive [2014/30/EU](#) of the European Parliament and of the Council of 26 February 2014 on the harmonisation of the laws of the Member States relating to electromagnetic compatibility (recast).
- Directive [2014/35/EU](#) of the European Parliament and of the Council of 26 February 2014 on the harmonisation of the laws of the Member States relating to the making available on the market of electrical equipment designed for use within certain voltage limits (recast).
- Directive [2011/65/EU](#) of the European Parliament and of the Council of 8 June 2011 on the restriction of the use of certain hazardous substances in electrical and electronic equipment and its amendments (Directive [2015/863/EU](#)).

Standards which Conformity is Declared:

- [EN 61326-1:2013](#) Electrical equipment for measurement, control and laboratory use - EMC requirements -- Part 1: General requirements (IEC 61326-1:2012)
- [EN 61010-1:2010](#) Safety requirements for electrical equipment for measurement, control and laboratory use -- Part 1: General requirements.

Barcelona, May the 1st 2017

A handwritten signature in black ink, appearing to read 'Diego Moreno', written over a horizontal line.

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IUL Inc.
IUL GmbH